

# Guide for People Data for TSI Link™ for Respiratory Protection



---

## Background

Ensuring clarity about what fields represent and how they are displayed in the system helps to ensure you import the data properly and understand what is shown in the Portal.

This guide provides links to an examples file, to an import template and provides a data dictionary. This can help you understand the People Data for TSI Link™ for Respiratory Protection.

---

## Links to Resources

[People Import Template](#)

[People Import Example](#)

[TSI Link for Respiratory Protection API documentation](#) (Login Required)

---

## Data Dictionary

Respirator Fields

Field Name	TSI Link field Name *	Required	Field Type	Example	Notes
lastName	Last Name	Yes	toText(64)	Storley	
firstName	First Name	Yes	toText(64)	Jill	
idNumber	Id Number	Yes	toText(64)	7878	Typically the employee number from the company, assigned by the company / customer
company	Company	No	toText(64)	Quest	
location	Location	No	toText(64)	Shoreview	
note	Note	No	toText(128)	Has some facial hair, take extra care to ensure proper fitting	text field that can be filled in with comments related to the person and their fit testing dynamics
custom1Label	<i>Custom Field 1 Label, name set by customer</i>	No	toText(64)	Department	Text field, null if not used
custom1Data	<i>Custom Field 1 value selected</i>	No	toText(64)	Marketing	Text field, null if not used
custom2Label	<i>Custom Field 2 Label, name set by customer</i>	No	toText(64)	Hazard Level	Text field, null if not used

Field Name	TSI Link field Name *	Required	Field Type	Example	Notes
custom2Data	<i>Custom Data 2 value selected</i>	No	toText(64)	Low	Text field, null if not used
custom3Label	<i>Custom Field 3 Label, name set by customer</i>	No	toText(64)	User Frequency	Text field, null if not used
custom3Data	<i>Custom Data 3 value selected</i>	No	toText(64)		Text field, null if not used
custom4Label	<i>Custom Field 4 Label, name set by customer</i>	No	toText(64)		Text field, null if not used
inactive	Active	No	toBoolean	FALSE	toBoolean: "yes", "true", 1="true", "no", "false", 0 = false
email	Email	No	toText(384)	david@tsi.com	
respiratorRequired	Respirator Required	No	toBoolean	TRUE	toBoolean: "yes", "true", 1="true", "no", "false", 0 = false
trainingDate	Last Respirator Training	No	toISODate	2024-06-26T18:40:43.355Z	must be an ISO date string "2025-01-01T08:00:00.000Z"
trainingExpiration	Respirator Training Expires	No	toISODate	2024-06-26T18:40:43.355Z	must be an ISO date string "2025-01-01T08:00:00.000Z"
isSupervisor	Is a Supervisor?	No	toBoolean	FALSE	toBoolean: "yes", "true", 1="true", "no", "false", 0 = false

Note \* Italics = not a specific field on the TSI Link website, so the use is described



Distributed by:  
 Kenelec Scientific Pty Ltd  
 1300 73 22 33  
 sales@kenelec.com.au  
 www.kenelec.com.au



**Knowledge Beyond Measure.**

TSI Incorporated – Visit our website [www.tsi.com](http://www.tsi.com) for more information.

USA Tel: +1 800 680 1220 India Tel: +91 80 67877200  
 UK Tel: +44 149 4 459200 China Tel: +86 10 8251 6588  
 France Tel: +33 4 91 11 87 64 Singapore Tel: +65 6595 6388  
 Germany Tel: +49 241 523030

TSI and the TSI logo are registered trademarks of TSI Incorporated in the United States and may be protected under other country's trademark registrations.

