# Setup Guide

System Setup and Configuration

**Vaisala viewLinc Monitoring System**

EN

DE

FR

ES

PT

ZH

JA

**VAISALA**

## PUBLISHED BY

Vaisala Oyj

Street address: Vanha Nurmijärventie 21, FI-01670 Vantaa, Finland

Mailing address: P.O. Box 26, FI-00421 Helsinki, Finland

Phone: +358 9 8949 1

Visit our Internet pages at [www.vaisala.com](http://www.vaisala.com).

No part of this manual may be reproduced, published or publicly displayed in any form or by any means, electronic or mechanical (including photocopying), nor may its contents be modified, translated, adapted, sold or disclosed to a third party without prior written permission of the copyright holder. Translated manuals and translated portions of multilingual documents are based on the original English versions. In ambiguous cases, the English versions are applicable, not the translations.

The contents of this manual are subject to change without prior notice.

Local rules and regulations may vary and they shall take precedence over the information contained in this manual. Vaisala makes no representations on this manual's compliance with the local rules and regulations applicable at any given time, and hereby disclaims any and all responsibilities related thereto.

This manual does not create any legally binding obligations for Vaisala towards customers or end users. All legally binding

obligations and agreements are included exclusively in the applicable supply contract or the General Conditions of Sale and General Conditions of Service of Vaisala.

This product contains software developed by Vaisala or third parties. Use of the software is governed by license terms and conditions included in the applicable supply contract or, in the absence of separate license terms and conditions, by the General License Conditions of Vaisala Group.

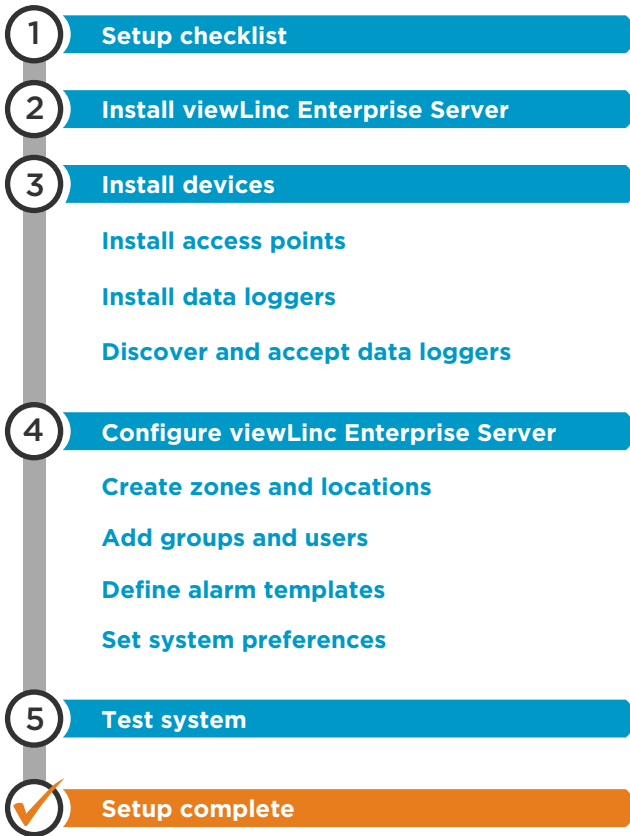
This product may contain open source software (OSS) components. In the event this product contains OSS components, then such OSS is governed by the terms and conditions of the applicable OSS licenses, and you are bound by the terms and conditions of such licenses in connection with your use and distribution of the OSS in this product. Applicable OSS licenses are included in the product itself or provided to you on any other applicable media, depending on each individual product and the product items delivered to you.

## Table of Contents

|                |    |
|----------------|----|
| English.....   | 5  |
| Deutsch.....   | 13 |
| Français.....  | 21 |
| Español.....   | 29 |
| Português..... | 37 |
| 中文.....        | 45 |
| 日本語.....       | 53 |

# Setup Roadmap

If you require assistance with the setup of your system, Vaisala Customer Care is ready to assist you, free of charge <sup>1)</sup> at [helpdesk@vaisala.com](mailto:helpdesk@vaisala.com).

1) Software product purchase entitles you to basic support, which is limited to hardware and software installation assistance. Comprehensive support is available through the purchase of a support plan.

# Setup Checklist

| viewLinc Enterprise Server |  |
|----------------------------|--|
| <input type="checkbox"/>   | I have a system administrator or IT support available to help me with network and server configuration.  |
| <input type="checkbox"/>   | A network server has been allocated to run viewLincEnterprise Server 24x7. I know its IP address and hostname, and it meets minimum installation requirements (see <i>viewLinc User Guide</i> or datasheet).   |
| <input type="checkbox"/>   | I know the address of the Network Time Protocol (NTP) server in the network.   |
| <input type="checkbox"/>   | My laptop or workstation has a supported web browser (Google Chrome™, Microsoft® Internet Explorer® 11, or Microsoft Edge™) and a PDF reader.  |
| <input type="checkbox"/>   | I know where the program files and viewLinc database should be stored. A backup system is in place to make sure the files are recoverable in case of server failure.   |
| <input type="checkbox"/>   | I know if the local network security policy requires me to use a trusted TLS certificate. If yes, I have located my company's signed certificate and key files, or have the information required to create a private certificate and key during installation (hostname, alias, number of years certificate will be valid for). |
| <input type="checkbox"/>   | I have the USB drive that contains the viewLinc setup files and license key.   |

| Devices                  |   |
|--------------------------|---|
| <input type="checkbox"/> | I have a plan that lists the serial number and installation location of each device.  |
| <input type="checkbox"/> | The plan includes information on how each device will be mounted (screws, magnets, etc).  |
| <input type="checkbox"/> | IP addresses have been allocated for my HMT330 series transmitters and data loggers' network connections (DHCP address reservations or static IPs). |
| <input type="checkbox"/> | Planned locations of wired devices are within 180 cm (6 feet) distance of power outlets.  |
| <input type="checkbox"/> | I know if drilling or door access is required for probes (for example, freezer installation).   |

| Network                  |  |
|--------------------------|--|
| <input type="checkbox"/> | Ethernet network drops with RJ45 jacks have been installed and tested at each location where a wired network connection is needed.     |
| <input type="checkbox"/> | A Power over Ethernet (PoE) injector or switch provides operating power at network drops where a PoE powered device will be installed. |

# Install viewLinc Enterprise Server Software

The viewLinc Enterprise Server software is delivered on a USB drive. Your license key is included with the drive.

For information on installation requirements, see *viewLinc User Guide*.

- ▶ 1. Insert the USB drive into your computer.
2. If you want to install viewLinc Enterprise Server on a remote server (not the local computer), copy the **viewLincInstall.exe** file from the USB drive to the target server.
3. Run **viewLincInstall.exe**. The installation wizard starts.
4. Follow the instructions on screen to complete the installation wizard. If an earlier version of the viewLinc software is already installed on the server, it will be upgraded to the new version. If the installation is an upgrade, some installation choices may be skipped or modified as necessary.
  - a. Select the installation language. The installation wizard will use the selected language, and it will be set as the system language for viewLinc Enterprise Server.
  - b. Enter your viewLinc License Key.
  - c. Select installation package type: **viewLinc Enterprise Server** or **Device Host** (additional servers for managing devices in a large installation).
  - d. Accept the license agreements.
  - e. Select destination folder for viewLinc software.
  - f. Select destination folder for viewLinc data.
  - g. Upload your TLS certificate and key, or have the installation wizard create a new certificate and key for you. If you are upgrading an old viewLinc version, you may also keep the existing certificate and key.
  - h. If the installation is an upgrade, you will be reminded to create a backup copy of your current viewLinc database. Always create a new backup before upgrading an existing viewLinc system.
  - i. Verify your selections and wait for the installation to complete.

viewLinc Enterprise Server software is now installed on your computer. The software will run in the background when the server is on, even when the user interface is not open.

# Install Devices

See device-specific documentation for detailed instructions.

If you are upgrading an existing system without adding new devices, you can skip this section.

Create a device installation plan to keep track of data logger serial numbers and installation locations.

## Install Access Points

Set up your **VaiNet access points** so that:

- They are able to connect to the Ethernet network
- They have the IP address or hostname of at least one NTP server they can access
- Their VaiNet channel number is set
- They have the IP address or hostname of viewLinc Enterprise Server
- They are in installation mode so that they can connect new VaiNet data loggers to the system

If you have new **WLAN access points**, configure their network and wireless settings.

When your wireless access points are set up, install them at their planned locations:

1. Mount the access point as described in its documentation.
2. Check that all cables are routed and secured appropriately.
3. Apply a location identification label to the access point according to your company policy.
4. Make sure the access point is turned on.

## Install Data Loggers and Transmitters

Set up your **data loggers** and **transmitters** so that they are ready for installation.

If you have **Vaisala vNet PoE Data Logger Interfaces** for your wired data loggers, follow vNet documentation to connect and configure them.

Mount the data loggers and transmitters at their planned locations:

- Mount the device as described in its documentation. Note that each device model has unique installation requirements.
- Apply a location identification label to the device according to your company policy.
- Make sure the device is turned on.

# Discover and Accept Devices

## New devices (3)

Data loggers must be accepted to the system before their data can be collected and used.

- ▶ 1. Open the viewLinc Enterprise Server user interface by double-clicking the desktop icon , or by entering the server's IP address and port in a web browser's address field.
2. Log in to viewLinc Enterprise Server as the default administrator user (user name **admin**, password **viewL!ncAdm1n**).
3. If this is the first time logging in as the admin user, a brief setup tour starts automatically. Complete the tour to familiarize yourself with system setup.
4. When viewLinc detects new devices, **New Devices** text appears at the top of the screen. Click it to open the **New Devices** window.
5. Compare the list of new devices to your device installation plan to verify that all your devices have been found. In the **Accept** column, select each device you want to add.
6. Select **Accept** and **Save** to register the selected devices with your viewLinc Enterprise Server.
7. Open the **Sites Manager** window and select the **Hosts and Devices** tab. On the **Hosts and Devices** tree, verify that all of your devices are now listed.

- If none of your devices are listed, check that network traffic to ports used by viewLinc (for example, 12600, 771, and 6767) is not blocked by a firewall.
- If some devices are missing, make sure they are configured, powered on, and connected to the network. Refer to the device-specific documentation for troubleshooting assistance.
- If the automatic discovery is not finding all of the devices, refer to the *viewLinc User Guide* for instructions on how to discover devices manually.

# Configure viewLinc Enterprise Server

After you verify that your devices are accepted and available, you are ready to configure the viewLinc Enterprise Server. Several interactive tours are available from the viewLinc desktop **Help > Tours** menu, to guide you through these configuration activities:

- Create Zones and Locations (watch a tour: Create a Zone/Create a Location)
- Add groups and users (watch a tour: Create a Group/Add a User)
- Add alarm templates (watch a tour: Create/Apply a Threshold Alarm Template; Create/Apply an Alarm Notification Template)
- Set system preferences (such as system language, measurement unit display, email/SMS settings)
- Optional setup tasks: assign permissions, set up schedules, create predefined comments

For more information about configuring viewLinc, refer to the instructions in the *viewLinc User Guide M211975EN*.

Download the **Planning Worksheet** (pdf) from **Overview > Getting Started** to define:

- Names for your monitored sites (Zones and Locations)
- Users, their assigned groups, rights and permissions
- Alarm thresholds required for different Locations
- Alarm notification delivery methods (email/SMS/command), and notification escalation paths

# Test the System

**System Test:** Complete a standard system test to ensure all components are installed and functioning as expected. Recommended for non-GxP installations. For testing instructions, see *viewLinc User Guide*.

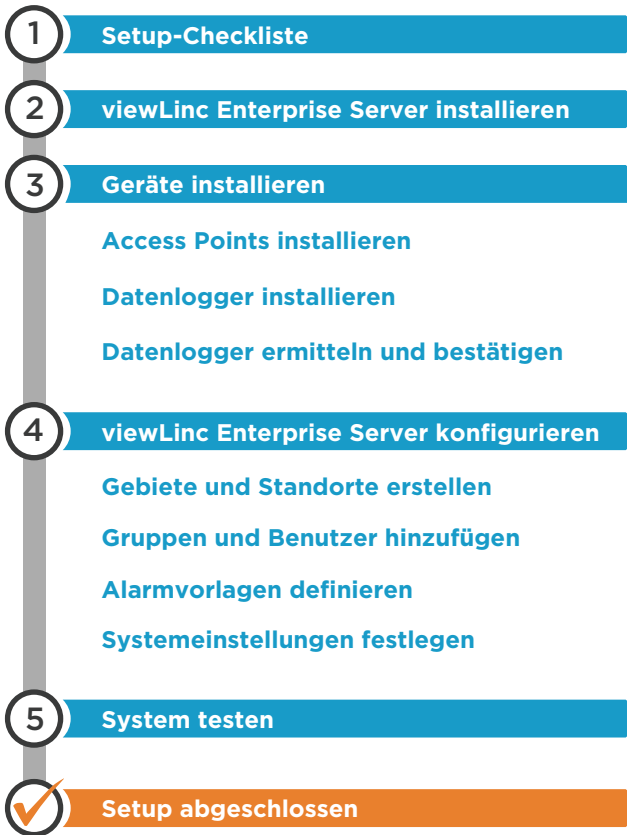
**IQOQ Protocol:** For controlled GxP environments, Vaisala provides a comprehensive **Installation Qualification and Operational Qualification Protocol** to ensure the system installation meets all your alarm monitoring requirements.

Contact Vaisala to obtain a viewLinc 5 protocol, available in 7 languages.

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Kenelec Scientific Pty Ltd