

# TrakPro™ v5 Data Analysis Software



## USB Quick Start Guide

This guide will help you to quickly begin using your TrakPro™ Data Analysis software for first time use with a personal computer and TSI® Equipment.

### 1. Software Installation

Download TrakPro™ v5 software from the TSI.com website in the [Service/Support](#) page. Once downloaded onto your PC, click the application file to begin installation. Follow the on-screen instructions to install the software.

### 2. Initial Start-Up of your TSI® Equipment

Plug the instrument into an AC power source or install the batteries. Press the **On/Off** button to turn the instrument on.



### 3. Communications Setup

After the instrument has completed its initial warm up, connect the USB cable from the instrument to your computer.

The “USB-A” connection to the computer, the “USB-B” or “USB mini-B” connection to the instrument.

### 4. Verify Communications

To verify successful Instrument-to-Computer communications:

Start the TrakPro™ software by double-clicking on the icon on the desktop.

For TrakPro™ v5 software (AM520 and AM520i only):

- Click on the **Devices** tab located on the left side of the application window. Verify that the Instrument and Serial Number information appear.

### 4. Verify Communications *(continued)*

For TrakPro™ v4.7.x software (all other TSI® products):

- Go to **Options | Software Configure...** menu. Choose the appropriate instrument for which you are connecting.
- Go to **Instrument Setup | Communications** menu. Verify that the Instrument and Serial Number information appear.

#### NOTICE

If the Instrument Model and Serial Number are not displayed, check driver installation by following the steps in the “TrakPro Data Analysis Software Installation Guide.”

### Further Learning

You are now ready to learn about the daily operations of your TSI® Instrument and the TrakPro™ software. Refer to the Operation and Service manual for your equipment and the TrakPro™ Data Analysis Software manual for detailed instructions on how to use the various functions of instrument and software.



Knowledge Beyond Measure.

TSI Incorporated – Visit our website [www.tsi.com](http://www.tsi.com) for more information.

USA	Tel: +1 800 680 1220	India	Tel: +91 80 67877200
UK	Tel: +44 149 4 459200	China	Tel: +86 10 8219 7688
France	Tel: +33 1 41 19 21 99	Singapore	Tel: +65 6595 6388
Germany	Tel: +49 241 523030		

P/N 6006952 Rev C

©2022 TSI Incorporated

Printed in U.S.A.



Distributed by:

Kenelec Scientific Pty Ltd  
1300 73 22 33  
[sales@kenelec.com.au](mailto:sales@kenelec.com.au)  
[www.kenelec.com.au](http://www.kenelec.com.au)