VAISALA

QUICK GUIDE

Vaisala Veriteq vLog

English
Français
中文
日本語

M211338EN-C

PUBLISHED BY

Vaisala Canada Inc
13775 Commerce Parkway
Richmond, BC V6V 2V4
Canada
1-866-861-3388 (or 604-273-6850)
veriteqsupport@vaisala.com

Visit our internet pages at www.vaisala.com

© 2011 Vaisala Canada Inc

No part of the manual may be reproduced in any form or by any means, electronic or mechanical (including photocopying), nor may its contents be communicated to a third party without prior written permission of the copyright holder.

The contents are subject to change without prior notice.

Please observe that this manual does not create any legally binding obligations for Vaisala towards the customer or end user. All legally binding commitments and agreements are included exclusively in the applicable supply contract or Conditions of Sale.

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English

Overview

This Guide is designed to get you started gathering data with vLog right away. It shows you how to install vLog software and points you to important information on setting up and using your Vaisala Veriteq data loggers.

Installation Requirements

Vaisala Veriteq vLog requires a PC with the following minimum configuration:

- 32-bit or 64-bit version of: Windows® XP, Windows Server® 2003, Windows Vista®, Windows Server® 2008, Windows® 7
- CD-ROM drive (if installing from a CD)

Quick Start Steps

Install vLog Software

- 1 Ensure you have administrator privileges.
- 2 Start Windows, log in as the administrator, and close all running applications. If you will be using

the USB Logger Cable, do not connect the cable until the installation is complete.

- 3 Insert the installation CD. The setup program starts automatically (if your system does not automatically run CD-ROMs, click **START** then **RUN**. If your CD-ROM drive is D, type d:\setup, then press **ENTER**.)
- 4 Select the language, then click **OK**. The vLog Setup wizard opens.
- 5 Click **NEXT**.
- 6 Accept the License Agreement, then click **NEXT**.
- 7 Enter your vLog security key, including spacers (for example, xxxx-xxxx-xxxx-xxxx-xxxx).
- 8 Select the destination folder, then click **NEXT**.
- 9 Accept or edit the location of the Program Folder (where vLog appears in Start Menu), then click **NEXT**.
- 10 If you want a desktop shortcut, select this option, then click **NEXT**.
- 11 Click **INSTALL**.
- 12 Click **FINISH**.

Connect Loggers to PCs

- Connect your logger to your PC using either a Serial Logger Cable, a USB Logger Cable, a vNet

or Digi device. For technical support, visit www.vaisala.com/lifescience-hitech, then select Service and Support.

Confirm the COM Port

Before you can communicate with your Vaisala Veriteq data logger, you need to know the number of the COM port that has been allocated for the data logger.

To check which COM port has been allocated, open Windows Device Manager:

- **Windows® XP:** Start | Settings | Control Panel | System | Hardware | Device Manager
- **Windows Server® 2003:** Start | Control Panel | System | Hardware | Device Manager
- **Windows Server® 2008/Windows Vista®:** Start | Control Panel | System | Device Manager
- **Windows® 7:** Start | Control Panel | Hardware and sound | Device Manager

Under Ports, note the COM port number allocated to the data logger.

Set up vLog to Communicate with the Data Logger

- 1 Select Start | Programs | Vaisala | Vaisala Veriteq vLog | Vaisala Veriteq vLog.

- 2 If you are running vLog VL, the first time you use vLog, the vLog Security Wizard opens. Click **NEXT**.

If you are using vLog SP, vLog starts automatically - go to step 9.

- 3 Enter the vLog security key you received with your package. Click **NEXT**.
- 4 Create a vLog administrator account. The user name and password you use for vLog must already exist on the PC.
- 5 Click **NEXT**. Specify the folder where you want to keep the audit trail file, either on the local computer or in a network location. You can modify this option later using **Tools | Administrator Options**.

The audit trail file is a log of operations performed in vLog by various users. If multiple PCs use the same loggers, they should use the same audit trail file.

- 6 Click **NEXT**.
- 7 Select the security level you want to apply to all vLog users (you can modify this option later).
- 8 Click **FINISH**. vLog starts.
- 9 From vLog, select **Tools | Options**, select the **General** tab, then choose the COM port allocated to the data logger. Click **OK**.

- 10 From vLog, select **Logger | Setup**. If communication with the data logger is established, the Logger Setup screen appears. You can use the Setup buttons on the right of the screen to configure the data logger, or, if you want to transfer data from the logger before making any changes, proceed to the next step.
- 11 To transfer data from the data logger, from vLog, select **Logger | Transfer**. Follow the instructions in the Transfer Wizard. After the transfer is complete, you will be asked to insert the .spl file into the active graph file. The .spl file extension refers to a logger file.
- 12 Select **OK** to insert the file into the graph, then select the channel data to insert, click **OK**.
- 13 Save using **File | Save**.

Technical Support

North America

Contact Vaisala Canada Inc, 8am-4pm PST Monday to Friday, at 1-866-861-3388 (or 604-273-6850) or email veriteqsupport@vaisala.com.

See also www.vaisala.com/lifescience-hitech.

For sales, pricing, quotations, or general information, please call 1-800-273-8374 (or 604-273-6850).

Vaisala Headquarters (Finland)

Vanha Nurmijärventie 21
01670 Vantaa
FINLAND
Industrial Instruments
Phone: +358 9 8949 2658
Fax: +358 9 8949 2295

Japan Service Center

42 Kagurazaka 6-Chome
Shinjuku-ku
Tokyo 162-0825
JAPAN

Phone: +81 3 3266 9611
Fax: +81 3 3266 9610

China Service Center

Floor 2, EAS Building
No. 21, Xiao Yun Road, Dongsanhuan Beilu
Chaoyang District
Beijing 100027
CHINA

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